

Alcogiene® NSAD70 RTU

Non-Sterile 70% Isopropyl Alcohol Disinfectant (Ready-to-Use)



Background

Alcogiene® NSAD70 Isopropyl Alcohol RTU disinfectants are designed to effectively eliminate microbial contamination in critical aseptic core manufacturing areas. Our high-quality products ensuring fast and thorough disinfection of surfaces and equipment. Whether you need to disinfect your entire cleanroom or target high-touch areas, our alcohol-based disinfectants provide the efficacy and convenience you need to achieve optimal hygiene levels. Alcogiene® NSAD70 Isopropyl Alcohol RTU disinfectants are available in both **Deionised Water** and **USP Grade Purified Water** formulations.

- ✓ 0.2 micron filtered and packaged into cleaned containers
- ✓ Filled and double bagged in a Grade C (ISO Class 7) cleanroom
- ✓ Fast evaporating and low residue
- ✓ Trigger sprayer bottles are fully assembled and ready-to-use
- ✓ Manufactured in a cGMP, HACCP and ISO 13485: 2016 Certified Facility

Variants Available

Variation

Isopropyl Alcohol 70%, Blended with DI

Isopropyl Alcohol 70%, Blended with USP Grade Purified Water

#ThinkHygieneThinkAlcogiene

Application Site

For use on walls, floors and hard surfaces.

Direction of Use

Thoroughly clean and removes any visible dirt or debris from the surfaces. For optimum results hold bottle upright and spray from a distance of 10cm to 20cm. Apply to surface to ensure complete coverage. Always close the nozzle after use. Allow surfaces to air-dry. No rinsing is required

Storage, Stability and Shelf Life

Store upright in original closed containers, away from sunlight and extremes of temperature. Store in a cool and well-ventilated place. Expiry date is printed on the product's label.

Features & Benefits

- 0.2 micron filtered and packaged into cleaned containers.
- Filled and double bagged in a Grade C (ISO Class 7) cleanroom.
- Fast evaporating and low residue.
- Manufactured in a GMP, ISO 9001 and ISO 13485 compliant facility.
- Trigger sprayer bottles are fully assembled and ready-to-use.
- Trigger sprayers offer stream delivery or coarse spray.

Efficacy Test

Efficacy Test	Contact time
EN 1276	5 min
EN 1650	15 min
EN 13697	15 mins
EN 16615	5 mins

Effective Against

- Aspergillus brasiliensis
- Candida albicans
- Enterococcus hirae
- Escherichia coli
- Pseudomonas aeruginosa
- Staphylococcus aureus



Packaging Size

Product Description

Alcogiene® NSAD70 Isopropyl Alcohol RTU
(Isopropyl Alcohol 70%, Blended with DI)

Size Quantity

Product Code

1 L 12/ctn ACG113IPA(70)DI-1.0TS

5 L 4/ctn ACG113IPA(70)DI-5.0P

200 L - ACG113IPA(70)DI-DRM

Alcogiene® NSAD70 Ethanol Denatured RTU
(Isopropyl Alcohol 70%, Blended with USP Grade Purified Water)

1 L 12/ctn ACG113IPA(70)UW-1.0TS

5 L 4/ctn ACG113IPA(70)UW-5.0P

200 L - ACG113IPA(70)UW-DRM

Distributed by :

Qualychem Solution PLT
Phone No.:
- 603 7831 0678 (Oversea customers)
- 1 700 818 105 (Malaysian ONLY)
Email:
- qualychemplt@gmail.com

QualyChem

XENIUM 仪翔科技

China

深圳市仪翔科技有限公司
联系方式: 0755-8952 2407